One step solution for cell culture and storage

# COS banker III-


COS Banker is a one-step solution for cell culture and storage. Suspend culture cells in COS Banker and store at -80 degrees for long term storage (Liquid nitrogen storage is not necessary). This is a serum-free solution with no animal products added.





### Features and Advantages

- Ready to Use (No Dilution Required !)
- Possible to store the Cells at -80°C in Deep Freezer
- Possible to retain Cells in high Viability after defrosting
- Does not contain Serum or Animal products\*
- Endotoxin-free tested / Bacteria free tested

\* "Cos Banker II" contains Human Serum Albumin (GMP Grade) as Stabilizer

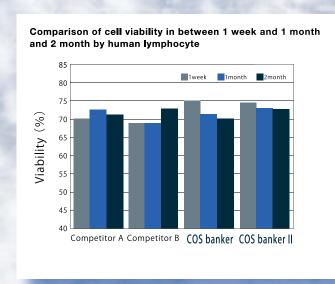

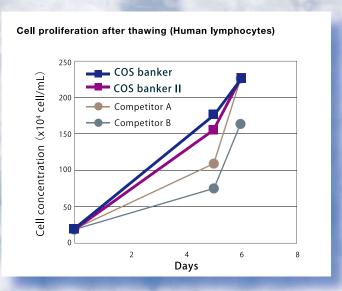

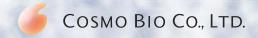

## One-step solution for cell culture and storage COS banker COS banker II



#### Instructions for preparation

- **1.** Collect culture cells (during logarithmic growth stage) and place  $5x10^5 1x10^7$  cells in a cryovial. Centrifuge for 5 minutes at 1,200 rpm.
- 2. Add 1 ml of Cos Banker to cell pellet and suspend. Freeze and store at -80°C.
- 3. If liquid nitrogen storage is desired, freeze cells at -80°C first before transfering to liquid nitrogen.
- 4. To thaw, defrost in a 37°C water bath quickly.

Note: It is ideal to use a logarithmic growth stage culture in order to increase the survival rate of cells when thawed.

#### **Tested Applications**

COS banker	■ Daudi (Human Burkitt lymphoma)	CTLL-2 (IL-2 depending mouse T-cells)
	■ HeLa S3 (Human cervical cancer)	■ NIH/3T3 (Mouse fatal cells)
	■ K562 (Human Chronic myelogenous leukemia)	■ PC-12 (Rat adrenal Pheochromocytoma)
	■ Raji (Human Burkitt lymphoma)	■ CHO-K1 (Chinese hamster ovarian cells)
	■ MOLT-4 (acute lymphoblastic leukemia: T-cells)	
COS banker II	■ Activated lymphocytes from Human Peripheral blood	

#### **Specifications**

Description	COS banker	COS banker II
Cat.No.	KOJ-COS-CFM01	KOJ-COS-CFM02
Size	120 mL	120 mL
Specification	Contains 10% DMSO without serum	Contains 10% DMSO and
		human serum albumin without FBS
Storage	2-8℃ in a cold dark place	2-8℃ in a cold dark place
Exp.	2 years after date of manufacture	2 years after date of manufacture

COSMO BIO CO., LTD.
TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU. TOKYO 135-0016, JAPAN

TEL: (81)3-5632-9617
FAX: (81)3-5632-9618
e-mail: export@cosmobio.co.jp
URL: www.cosmobio.com



10189