



Collagen coating solution

Catalog No. SCO

May 18, 2021

Catalog No.	Product Name	Size	Storage Conditions	Expiration date
SCO	Collagen coating solution (Type I collagen, Bovine)	100 mL	4°C	- Written on the bottle

Materials required

- ◇ Culture dish
- ◇ Collagen coating solution(#SCO)
- ◇ PBS (-) or HBSS (-)

Collagen Coating.

1. Put the collagen coating solution (PMC, Code No.: SCO) in a culture vessel to be used at a degree of being covered, leave the resulting material to stand for one hour or more (overnight for the glass bottom), and then remove the solution.
2. Rinse the vessel with PBS twice, and then put PBS therein, and leave the resultant material to stand for one or more hours, and then use the resultant material. If the material is not used immediately, fill the vessel with PBS, which can be kept for several days at room temperature.



COSMO BIO Co., LTD.

【JAPAN】
TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU. TOKYO 135-0016, JAPAN
Phone: +81-3-5632-9610
FAX: +81-3-5632-9619
URL: <https://www.cosmobio.co.jp/>



COSMO BIO USA

【Outside Japan】
2792 Loker Ave West, Suite 101
Carlsbad, CA 92010, USA
email: info@cosmobiousa.com
Phone/FAX: (+1) 760-431-4600
URL: www.cosmobiousa.com



コスモ・バイオ株式会社

COSMO BIO CO., LTD.

For research use only. Not for clinical diagnosis.

コラーゲンコート用溶液 (Collagen coat solution, Code No.SCO)

2020年5月18日訂正

※本マニュアルをご精読の上、研究目的にのみご使用ください。

Catalog No.	Product Name	Size	Storage Conditions	Expiration date
SCO	コラーゲンコート用溶液 (ウシ由来 I 型コラーゲン)	100 mL	4°C	-ボトルに記載

【事前に準備するもの】

- ・ご使用になる培養容器
- ・コラーゲンコート用溶液(製品コード:SCO)
- ・PBS(-)または HBSS(-)

コラーゲンコート方法

※コラーゲンコート用溶液は薄めずにそのままお使いいただけます。

1. 使用する培養容器にコラーゲンコート用溶液(製品コード SCO)が覆われる程度加え(24 ウェルプレートの場合 300~500 μ L)、1 時間以上(ガラス底の場合は一晩)静置後に溶液を除去します。
2. PBS で 2 回洗浄後、PBS を加えて 1 時間以上静置してからご使用してください。ただちに使用しない場合は PBS を満たして室温で数日間保管可能です。



コスモ・バイオ株式会社
COSMO BIO Co., LTD.

〒135-0016 東京都江東区東陽 2-2-20 東陽駅前ビル
URL : <http://www.cosmobio.co.jp/>

● 営業部（お問い合わせ）

TEL : (03) 5632-9610 FAX : (03) 5632-9619
TEL : (03) 5632-9620

● 札幌事業部（技術的なお問い合わせ）

TEL : (0134) 61-2301 FAX : (0134) 61-2295
E-mail : primarycell@cosmobio.co.jp