FITC-conjugated anti-Collagen IV α 5(IV) chain, Human (Mono)

FITC-Anti Collagen IV α 5(IV) Chain, Human (Mono) + Texas Red-Anti Collagen IV α 2(IV) Chain, Human (Mono)

Cat. No. SGE-CFT453

www.cosmobio.co.jp

Code: SG-CFT453

Quantity: 500 μ l

Appearance: Solution. Monoclonal antibody was purified by affinity chromatography.
The monoclonal antibody against α 5(IV) was conjugated with FITC.
0.1% NaN_3 is added for preservation.

Clone names: H53 (rat IgG2a/kappa)

Specificity of antibodies: H53 is specific to imperfection III of human α 5(IV).
Not reactive with human α 1, α 2, α 3, α 4 and α 6.
The amino acid sequence of the epitope is IDVEF. (Reference 1)

Conjugate : FITC


Preparation of antibodies: Monoclonal antibody was prepared by the rat lymph node method developed by Shigei Medical Research Institute with synthetic peptide of type IV collagen as immunogen.

Use: Staining of human cryostat sections by direct immunofluorescence (No acid treatment is necessary.)

Storage: Store in dark, 2-4°C, stable 2 year under this condition.

Reference: 1) Kagawa *et al.* (1997) *Nephrol. Dial. Transplant.* **12**: 1238-1241

Producer: Shigei Medical Research Institute

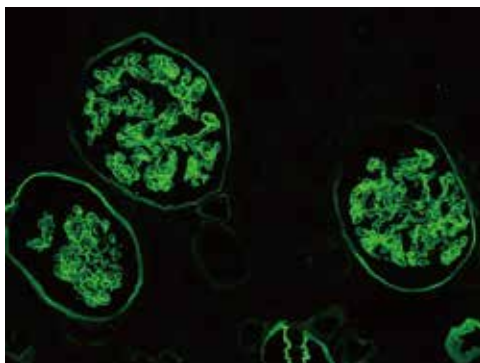The chromatic figure of normal human kidney stained with the product

As the procedure is direct-staining without acid treatment and possible to obtain the staining result easily and rapidly. Applicable to staining for basement membrane of glomerular, Bowman's capsular and partial renal tubular.

For research use only, Not for diagnostic use.
Please read this manual thoroughly before use.

Manufacturer

Shigei Medical Research Institute

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COSMO BIO CO., LTD.