


Instructions for stool collection for DNA stabilization (Spoon type)

☆ Please check the contents of kit.

Collection container (containing liquid): 1 tube
Stool collection sheet: 1 sheet
(This sheet can be flushed after use.)
Instructions you are reading now: 1 sheet

How to collect stool


- ① Read instructions and directions for use printed on the stool collection sheet carefully before use.
- ② Move the bowels on the stool collection sheet.
- ③—1 Hold the spoon of the collection container downward to prevent leakage of the liquid in the tube.
—2 Pull out a spoon part by turning a brown cap.
—3 Please pay attention to prevent spilling the liquid.

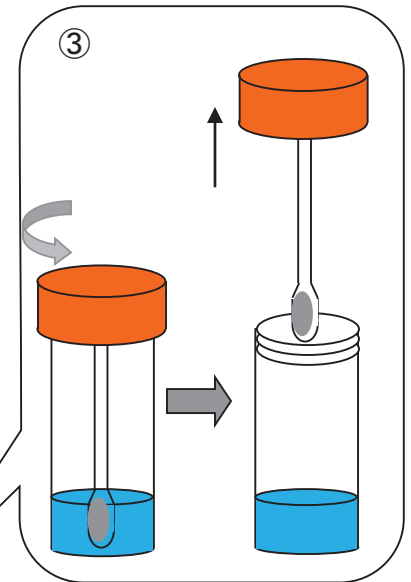

☆ Notice !!

In case of contact with the liquid (guanidine solution) in the container by mistake, wash in water immediately although the liquid causes no harm instantly. Please pay attention to children to prevent drinking it by mistake.

- ④ Collect stool by scooping it from several different places.

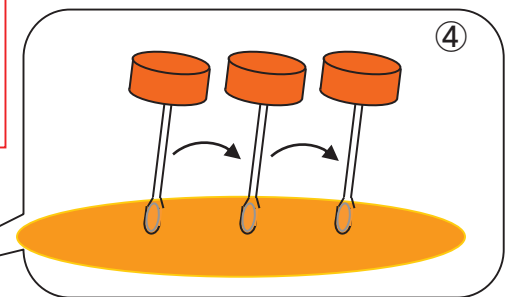


☆ Notice !!


(Adzuki bean grain size (0.5~0.7mm) of stool after several stabbings is sufficient amount.)

(Be careful to avoid protruding stool sample from the container due to too much stool.)

- ⑤—1 Return the spoon part into the collection tube with a liquid.
—2 Turn and tighten a brown cap to the end. Please make sure proper closure.
—3 Shake the container 5 or 6 times to mix the contents.

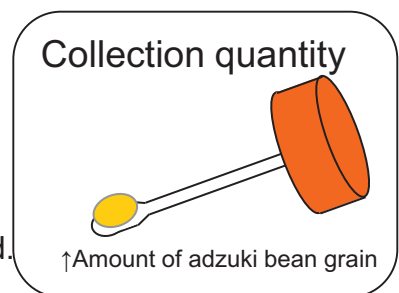

- ⑥ Please store the container with the sample in the supplied plastic bag at room temperature until the submission day.

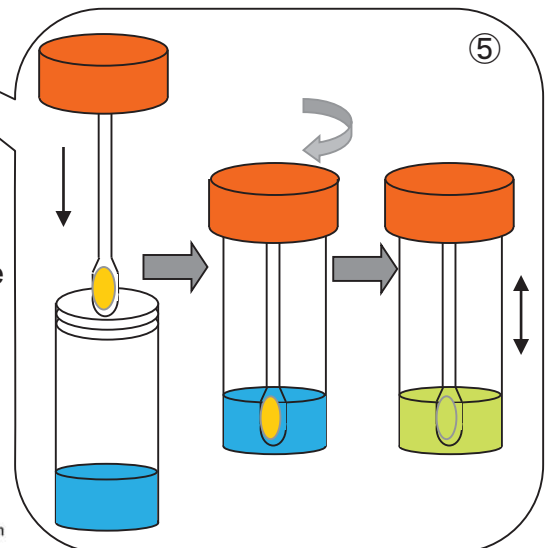

COSMO BIO CO., LTD.
Inspiration for Life Science

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME, KOTO-KU, TOKYO 135-0016 JAPAN
TEL: (81)3-5632-9617 / FAX: (81)3-5632-9618 / e-mail: export@cosmobio.co.jp / URL: www.cosmobio.com